Isotopx

Technical Brief T20812

Measurement of 2ng Nd (NdO⁺) using Xact Faraday collectors on a Phoenix TIMS equipped with $1e^{11} \Omega$ and $1T \Omega$ resistors.

Zenon Palacz, Davydd Wanless, Ian Turner & Tony Jones. Isotopx Ltd. UK

Introduction

We have shown that very small ion beams of Nd⁺ (5 to 50mv) can be measured accurately and precisely using the **Phoenix TIMS** equipped with either $1e^{11} \Omega$ or $1T \Omega$ resistors¹. For small sample sizes measurement of Nd as Nd⁺ is not the best approach since the ionization efficiency is not sufficiently high.

On the **Phoenix,** ion/atom efficiency for Nd⁺ is typically 2% to 2.5% across all sample sizes. It has been shown that measuring Neodymium as an oxide species provides higher ionisation efficiency than the metal and allows the routine measurement of Nd at the sub 10ng level².

In this work we have analysed 2ng aliquots of Nd as NdO using both $1e^{11} \Omega$ and 1T resistors to see if the lower noise characteristics of the 1T resistors provides an improvement in precision and accuracy. We have used the loading method of Harvey and Baxter² (2009) where the source of oxygen is Ta_2O_5 loaded with the sample. Much of the sensitivity produced by this method can be attributed to the loading protocol. Note that we do not believe we have perfected this method but this work is intended to show a comparison in the performance of the two types of resistor for small ion beam measurements, rather than the ultimate sensitivity and detection limit.

Sample Preparation

2ng of Jnd-i Nd standard was loaded onto single degassed rhenium filaments together with a Ta_2O_5 solution. The solutions were dried down at 0.5A, then heated slowly to a dull red colour for 2 seconds. Essentially the samples were loaded in the same way as for Strontium isotope ratio analysis.

Mass Spectrometry.

Filaments were loaded into a **Phoenix TIMS** spectrometer. Following pump-down the source vacuum was $1e^{-8}$ mbar and the analyser vacuum was $1e^{-9}$ mbar.

The filament was heated over a period of 40 minutes and an ion beam of 200 mv ($2e^{-12}$ amps) was optimized. Isotope ratios were measured in the dynamic mode in blocks of 20 cycles. Each cycle comprised 3 mass positions of 10 second integration. A delay of 3 seconds was made between mass jumps to accommodate the 3 second time decay to reach baseline for the 1T resistors¹. Typically 10 blocks of data were taken. It was found that the ion beam could last for much longer, however, the mass fractionation became too large for adequate correction. Baselines and ion focussing occurred at the start of each block of data. Baselines were measured at +/-0.5 amu for 30 seconds either side of each of the 3 mass positions. The same method was used for both $1e^{11}\Omega$ and 1T resistors.

Oxygen corrections were performed on-line using the oxygen isotope ratios of Nier 1950. It is known that this may not be totally accurate (Harvey and Baxter 2009²).

Results

Sensitivity and Mass Fractionation

Data for $1e^{11} \Omega$ and 1T resistors are shown in Table 1.

During analysis the ion signal remained fairly constant over the course of 3 to 3.5 hours, but towards the end of the measurement rapid increases and decreases in intensity occur (Figure 1).

The ion/atom efficiency obtained during analysis is typically 10%, though efficiencies of up to 17% are possible if the sample is taken to exhaustion.

The mass fractionation increases steadily over the course of a typical 3-4 hour analysis, but increases rapidly towards the end of the measurement (Figure 2). The exponential fractionation corrected ratio remains stable during the analysis (Figure 3). However more extreme mass fractionations are possible resulting in a non-exponential fractionation correction, typically when ¹⁴⁶Nd/¹⁴⁴Nd went above 0.7245. These data were not used in the compilation. Modifications in loading method may reduce the rate of mass fractionation.

Isotope ratios

Isotope ratio data for $1e^{11} \Omega$ and 1T resistors on 2ng aliquots of Nd are shown in Table 1. Comparative data for ¹⁴³Nd/¹⁴⁴Nd and ¹⁴⁵Nd/¹⁴⁴Nd are shown in Figures 4 and 5 respectively. Taken as a whole, there is no significant difference between the two data sets. This supports previous work¹ which showed no significant difference in the performance of 1T and $1e^{11} \Omega$ resistors for ion signals >50mV.

For the ¹⁴³Nd/¹⁴⁴Nd, internal precision is typically 5-7ppm 1RSE, external precision is in the range 15 to 20ppm 1RSD.

The 1T data are closer to the expected ratio of 0.512104 which is obtained for large loads (>100ng) on triple filaments measured as Nd⁺. This difference is likely to be related to the choice of oxygen isotope ratio.

The data are about two times less precise than NdO⁺ data reported by Harvey and Baxter 2009², though they used 4ng of Nd in their standard data set.

Conclusions

- Analysis of Nd as NdO+ is at least 5 times more sensitive than analysing as Nd+.
- 2ng loads of Nd measured as NdO+ can be analysed to internal precision of 7ppm 1rse and external precision of about 15ppm 1RSD.
- There is no difference between the performance of 1T and 1e¹¹ Ω resistors for Nd measured at these ion intensities.

Technical Brief T20812



2ng NdO 1T resistor data											
	¹⁴² Nd/ ¹⁴⁴ Nd	%1se	¹⁴³ Nd/ ¹⁴⁴ Nd	%1se	¹⁴⁵ Nd/ ¹⁴⁴ Nd	%1se	¹⁴⁸ Nd/ ¹⁴⁴ Nd	%1se	¹⁴² Nd ¹⁶ O amps	Minutes	ion/atom efficiency
1	1.141818	0.0008	0.512085	0.0006	0.348413	0.0006	0.241589	0.0012	3.3E-12	152	9.0%
2	1.141821	0.0008	0.512099	0.0008	0.348420	0.0009	0.241589	0.0015	2.9E-12	187	9.5%
3	1.141816	0.0007	0.512103	0.0007	0.348419	0.0007	0.241594	0.0010	3.1E-12	206	11.4%
4	1.141764	0.0010	0.512085	0.0007	0.348409	0.0006	0.241603	0.0013	3.1E-12	207	11.5%
5	1.141838	0.0010	0.512080	0.0007	0.348406	0.0006	0.241590	0.0015	3.1E-12	229	12.6%
6	1.141776	0.0010	0.512093	0.0007	0.348428	0.0006	0.241598	0.0015	3.1E-12	206	11.5%
7	1.141789	0.0009	0.512091	0.0007	0.348423	0.0006	0.241597	0.0014	3.2E-12	206	11.5%
8	1.141773	0.0011	0.512078	0.0007	0.348412	0.0007	0.241607	0.0016	3.1E-12	187	10.1%
9	1.141818	0.0009	0.512097	0.0007	0.348419	0.0006	0.241593	0.0013	3.2E-12	206	11.6%
10	1.141762	0.0014	0.512084	0.0009	0.348417	0.0008	0.241587	0.0014	3.2E-12	103	5.8%
11	1.141838	0.0007	0.512109	0.0006	0.348410	0.0006	0.241598	0.0011	3.0E-12	204	11.0%
Mean	1.141801	0.0009	0.512091	0.0007	0.348416	0.0007	0.241595	0.0013	3.1E-12	190	10.5%
1SD	0.000029	0.0002	0.000010	0.0001	0.000006	0.0001	0.000006	0.0002	1.1E-13	35	2%
1RSD	0.0025%		0.0019%		0.0018%		0.0026%				

2ng NdO 1e ¹¹ ohm resistor											
									¹⁴² Nd ¹⁶ O		ion/atom
	¹⁴² Nd/ ¹⁴⁴ Nd	%1se	¹⁴³ Nd/ ¹⁴⁴ Nd	%1se	¹⁴⁵ Nd/ ¹⁴⁴ Nd	%1se	¹⁴⁸ Nd/ ¹⁴⁴ Nd	%1se	amps	Minutes	efficiency
1	1.141797	0.0007	0.512084	0.0007	0.348407	0.0007	0.241579	0.0013	3.1E-12	184	10.3%
2	1.141790	0.0007	0.512072	0.0007	0.348400	0.0007	0.241585	0.0013	3.2E-12	179	10.1%
3	1.141825	0.0005	0.512079	0.0005	0.348416	0.0005	0.241598	0.0013	3.2E-12	162	9.1%
4	1.141812	0.0005	0.512076	0.0005	0.348408	0.0005	0.241592	0.0012	3.2E-12	181	10.1%
5	1.141838	0.0006	0.512084	0.0006	0.348397	0.0006	0.241598	0.0012	3.3E-12	181	10.5%
6	1.141774	0.0005	0.512062	0.0005	0.348404	0.0005	0.241601	0.0012	3.2E-12	182	10.5%
7	1.141817	0.0006	0.512081	0.0006	0.348400	0.0006	0.241579	0.0014	3.1E-12	183	10.0%
8	1.141804	0.0007	0.512068	0.0007	0.348401	0.0007	0.241577	0.0013	3.3E-12	174	10.1%
9	1.141786	0.0006	0.512076	0.0006	0.348387	0.0006	0.241585	0.0014	3.1E-12	183	9.9%
10	1.141853	0.0006	0.512084	0.0006	0.348406	0.0006	0.241574	0.0012	3.3E-12	179	10.4%
Mean	1.141810	0.0006	0.512077	0.0006	0.348403	0.0006	0.241587	0.0013	3.2E-12	179	10.1%
1SD	0.000025	0.0001	0.000007	0.0001	0.000008	0.0001	0.000010	0.0001	7.4E-14	6	0.4%
1RSD	0.0022%		0.0015%		0.0022%		0.0041%				

Table 1. Isotope ratios determined on 2ng NdO using 1e¹¹ ohm and 1T resistors.

Isotopx

Technical Brief T20812



Figure 1. Typical ion beam behaviour during the analysis of 2ng NdO



Figure 2. Mass Fractionation behaviour with time during analysis of 2ng NdO



Figure 3. Relationship between fractionation corrected ¹⁴³Nd/¹⁴⁴Nd and the normalizing ratio ¹⁴⁶Nd/¹⁴⁴Nd during analysis.

Isotopx

Technical Brief T20812



Figure 4. Comparative behaviour between 1T and 1e¹¹ ohm resistors for ¹⁴³Nd/¹⁴⁴Nd



Figure 5. Comparative behaviour between 1T and 1e¹¹ ohm resistors for ¹⁴⁵Nd/¹⁴⁴Nd

References

- Multidynamic measurement of small Neodymium (Nd⁺) ion signals on Faradays using X act amplifiers with 1e¹¹Ω and 1e¹²Ω (1T) resistors. Zenon Palacz, Davydd Wanless, Ian Turner & Tony Jones. Isotopx Technical Brief T10712, 2012. <u>Click to download.</u>
- An improved method for TIMS high precision neodymium isotope analysis of very small aliquots (1–10 ng). Harvey, J., Baxter, E.F. Chemical Geology Vol. 258 pp. 251–257, 2009. <u>Link to open access pdf</u>
- 3. A re-determination of the relative abundances of the isotopes of carbon, nitrogen, oxygen, argon and potassium. Nier, A.O., Physical Review, Vol. 77, Issue 6, pp. 789-793, 1950. Link to pdf, purchase required.