

How can SIRIX be used to improve CO₂ measurement?

Introduction

The objective of CO₂ isotopologue studies is to accurately and precisely measure small differences ($\Delta 47$) in the abundance of m/z 47 (¹³C¹⁸O¹⁶O) relative to m/z 44 ¹²C¹⁶O¹⁶O. Small changes in $\Delta 47$, are related to changes in temperature of formation of carbonate rocks¹. The relationship between temperature and $\Delta 47$ can be defined by the equation¹:

$$\Delta 47 = 0.0592 \times 10^6 \times T^{-2} - 0.02$$

Where T is the Temperature in Kelvin and $\Delta 47$ is in per mil (1/1000). This is depicted in Figure 1 below.

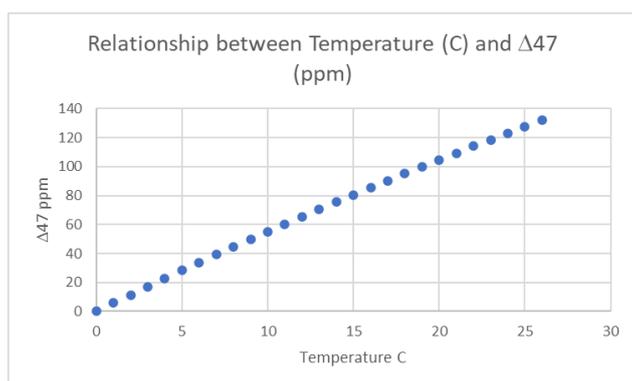


Figure 1. Relationship between $\Delta 47$ and temperature in carbonates. A resolution of 5 ppm will resolve a temperature difference of 1°C.

The isotope ratio 47/44 is approximately 4.5e-5. The mass spectrometric challenge to achieve these precisions include:

- A highly stable ion source
- High sensitivity from the ion source
- High resolution ion optics
- Flexible collector configuration to allow continuous baseline measurement
- No baseline aberrations from ion collisions with the flight tube
- Low noise, high dynamic range Faraday amplifier



Calcium carbonate tufa towers at Mono Lake in California

SIRIX large radius stable isotope ratio mass spectrometer

The SIRIX mass spectrometer design is essentially a Phoenix thermal ionization mass spectrometer, with the Thermal Ionisation source removed and replaced with a high efficiency electron bombardment gas source.


The SIRIX ion source is constructed of high purity stainless steel. It is pumped with a 300 l/s turbomolecular pump backed with an oil-free scroll pump. Ions are accelerated through a potential of 10 KV, through an exit slit of 0.3 mm width. Ions are focused in the X-Y-Z direction.

The large electromagnet produces a mass dispersion of 610 mm, this was designed for TIMS applications, but is well suited for CO₂ isotopologue studies as the isotope peaks are well separated at the collector, thereby reducing peak tailing. The base mass resolution of the SIRIX is 600 (10% valley) which is higher than most other gas source instruments. The high resolution ensures that hydrocarbons cannot interfere with peak centres.

An example of the peak shape and collector coincidence is shown in Figure 2. All collectors have 1mm entrance slits resulting in the same peak widths across the collector. Flat topped peaks are generated because the ion beam width is 0.3 mm.

Measurement of clumped CO₂ isotopologues using SIRIX stable isotope ratio mass spectrometer

AN22_02 / David Rousell

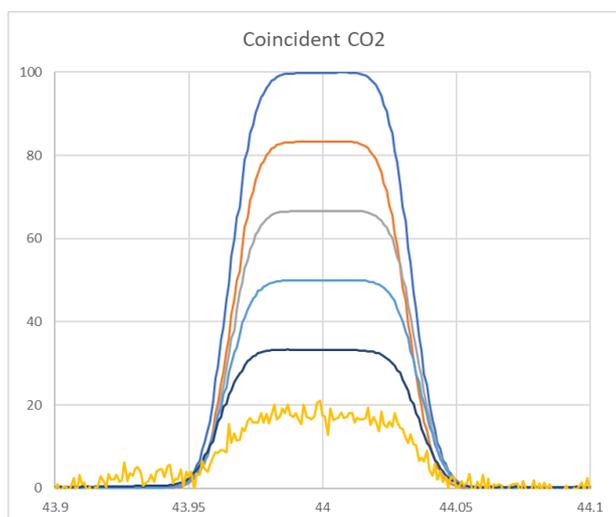Figure 2. Peak coincidence of CO₂ isotopologues using SIRIX

Flexible ion collection

A feature of the SIRIX is that all nine Faraday collectors can be moved independently. There is no fixed collector spacing. Thus no compromise in peak positioning and each isotope system is ensured of perfect peak coincidence.

In the context of CO₂ isotopologue measurements, not only can the isotopologues of CO₂ be measured in STATIC multicollection mode, but at the same time the half-mass baselines can be measured between the minor isotope peaks. This is shown in Table 1 below.

Collector	Mass	Typical intensity (amps)
H4	49	2.10E-14
H3	48.5	
H2	48	6.02E-13
H1	47.5	
Ax	47	6.38E-12
L2	46.5	
L3	46	5.60E-10
L4	45	1.54E-09
L5	44	1.40E-07

Table 1. Collector positions for CO₂ with an indication of the large range of ion intensities.

Baseline (peak tailing) measurement

The importance of accurate baseline measurement is demonstrated in Figure 3, which shows the high mass peak tailing from the *m/z* 44 peak beneath the minor isotopes 47, 48 and 49. The 44 signal is 1.4e-7 amps. The 47 peak is off-scale at 6.3e-12 amps. The peak tail at mass 47 is 5e-15 amps, so without a tail correction the contribution of the peak tail under mass 47 is approximately 0.1%.

At mass 48 the uncorrected peak tail is approximately 1% of the 48 peak. There is no evidence of any other baseline artefacts, for example from collisions with the flight tube. These reflections would produce larger errors in isotope ratios with increased ion intensity and would prevent an improvement in precision with increasing ion intensity. We believe the lack of reflections is due to internal baffling and the wide flight tube.

The amount of peak tail can be seen relative to the *m/z* 49 peak in Figure 2. The 49 peak is 0.14 ppm of mass 44. Therefore, at mass 47 it is 0.07 ppm relative to 44, at mass 48 it is 0.03 ppm and <0.01 ppm at mass 49. All the peaks are clearly resolved from the peak tail.

Table 1 shows that for this application the peak tail at mass 47 can be exactly quantified by measuring *m/z* 46.5 and 47.5. By taking the average of the two intensities a precise baseline correction can be made. For mass 48 the baseline is measured at 47.5 and 48.5. For 49 a baseline at 48.5 is used. Baseline subtraction can be made for each integration. This is important because during analysis the ion signal is constantly reducing, and baselines taken at the start of the measurement are not applicable during the measurement.

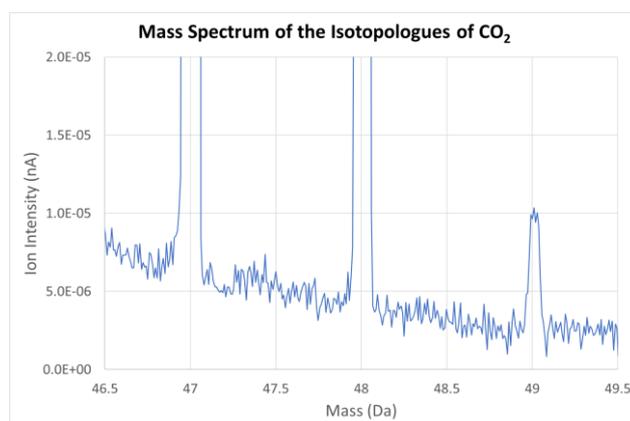


Figure 3. The high mass peak tail from mass 44 beneath 47, 48 and 49. No electrostatic filters are needed to remove the peak tail as the baseline can be corrected for with each ratio.

ATONA amplifier technology

The ATONA™ amplifier technology has been described elsewhere. In the example of isotopologue measurements, the ATONA is particularly well suited. The extremely large ion intensities of the major isotopes (Table 1) can be measured using the same detectors as the minor isotopes without range changing or calibration. Features of the ATONA include:

- Extremely rapid response; maximum signal to baseline in less than 300 ms
- Dynamic range of **ten** orders of magnitude
- Gain stability of <2 ppm 2RSD over 24 hours
- Noise stability of ~1e-18 amps over an hour.

CO₂ measurements

Figure 4 Shows typical ion beam behavior over the course of 100 minutes. Each data point is a 1 second integration. A total of 6000 integrations were made, in 10 blocks of 600. The ion beam was peak centered at the start of each block. The ion signal of the 44 CO₂ starts at just over 1000 volts (relative to a 1e11 ohm resistor) or 1e-8 amps. It decays slowly over the period by about 10%. No ion focusing was done during the acquisition. The mass 47 isotope was 0.4V, or 4e-12 amps. This is too high for a 1e13 ohm resistor but could be measured with a 1e12 ohm resistor.

The 45/44 isotope ratio shows some fine structure during the acquisition, and this is also seen within the 47/44 data set. We believe this is due to mass fractionation effects within the ion source.

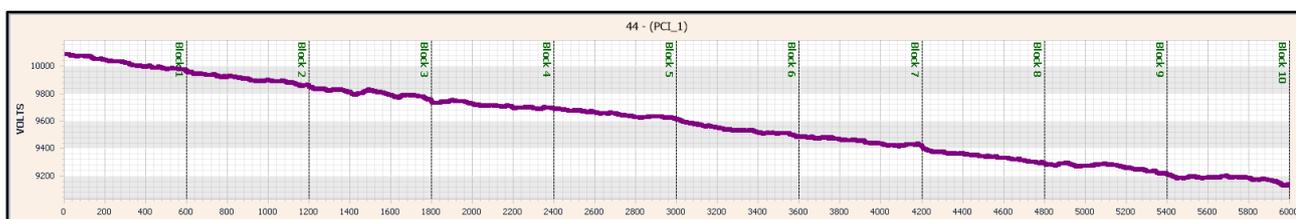


Figure 4: m/z 44 intensity as a function of time over 100 minutes

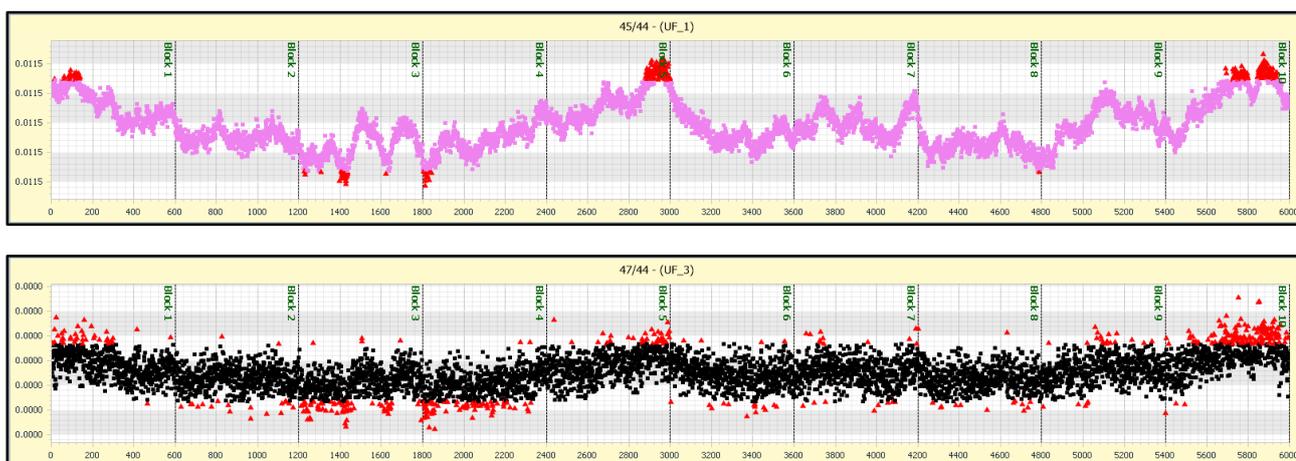


Figure 5: 45/44 and 47/44 ratios over the same time period.

Measurement of clumped CO₂ isotopologues using SIRIX stable isotope ratio mass spectrometer

AN22_02 / David Rousell



The final results are shown in Table 2 below.

	Mean	1se (ppm)
45/44	1.1487E-02	0.77
46/44	4.0674E-03	1.49
47/44	4.5230E-05	3.2
48/44	4.2066E-06	13.2
49/44	7.0171E-08	544

Table 2. Final data with mean values and precision values

Results

The results can be further broken down into individual 10 minute, 600 cycle integrations. This is shown in Table 3. Individual data blocks produce 47/44 data of sub-10 ppm 1rse, while the major isotopes have precisions of less than 2 ppm 1rse. The reproducibility expressed as 1RSD is not as good as this. For example, the reproducibility of 45/44 is 54 ppm, for 46/44 it is 107 ppm and for 47/44 it is 168 ppm.

The increase is entirely related to the mass difference between the isotopes and not the absolute abundance of the isotope. Despite the 47 isotope being two orders of magnitude smaller than the 46 isotope, the reproducibility of the data is controlled by mass fractionation. This is further illustrated in Figure 5, where a strong correlation is shown between 47/44 and 45/44 and follows a trajectory of mass fractionation with the loss of the lighter 44 isotope. The corresponding data summary is shown below.

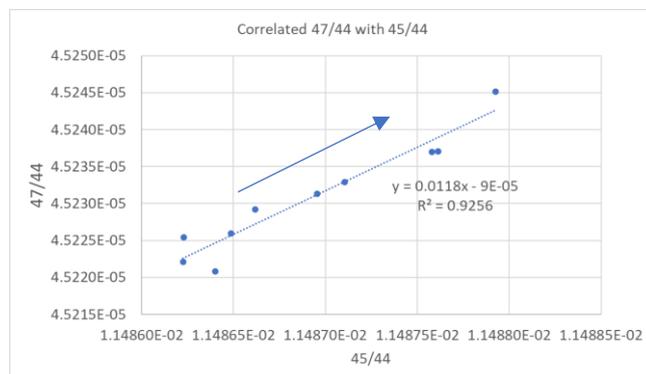


Figure 5. Correlation between 47/44 and 45/44 over 2 hours of data collection. Early data are at the lowest end of the trajectory with the arrow showing increasing time.

Summary

SIRIX can measure all the isotopes of CO₂ with high precision and accuracy. Large ion signals in excess of 1000 volts relative to a 1e11 ohm resistor (>1e-7 amps) can be measured with ATONA. Peak tailing from mass 44 is not a significant source of error. 47/44 precisions of ~3 ppm 1rse can be generated. Higher precision is limited to accurate correction of mass fractionation produced in the ion source.

Reference

1. P. Ghosh, J. Adkins, H. Affek, B. Balta, W. Guo, E. Schauble, E.D.Schrag. 13C-18O bonds in carbonate minerals: A new kind of palaeothermometer. *Geochemica et Cosmochimica Acta* 2006. 70. 1439.

TIME (mins)	Block	Intensity in Vots relative to 1e11 ohm resistor						Isotope ratios									
		44	45	46	47	48	49	45/44	ppm 1se	46/44	ppm 1se	47/44	ppm 1se	48/44	ppm 1se	49/44	ppm 1se
10	1	10,027	115.2	40.8	0.454	0.042	0.00070	1.1488E-02	1.8	4.0680E-03	3.5	4.5237E-05	8.8	4.2076E-06	42	7.0221E-08	1741
20	2	9,910	113.8	40.3	0.448	0.042	0.00069	1.1486E-02	0.8	4.0672E-03	1.5	4.5226E-05	7.6	4.2057E-06	43	7.0011E-08	1712
30	3	9,806	112.6	39.9	0.443	0.041	0.00069	1.1486E-02	1.8	4.0670E-03	3.5	4.5222E-05	9.1	4.2053E-06	41	6.9918E-08	1721
40	4	9,717	111.6	39.5	0.439	0.041	0.00068	1.1486E-02	1.6	4.0671E-03	2.9	4.5221E-05	9.0	4.2056E-06	42	6.9896E-08	1562
50	5	9,652	110.9	39.3	0.437	0.041	0.00068	1.1488E-02	2.1	4.0679E-03	3.8	4.5237E-05	9.1	4.2075E-06	40	7.0226E-08	1746
60	6	9,541	109.6	38.8	0.432	0.040	0.00067	1.1487E-02	1.3	4.0673E-03	2.5	4.5229E-05	8.7	4.2065E-06	39	7.0072E-08	1711
70	7	9,453	108.6	38.4	0.428	0.040	0.00066	1.1487E-02	1.3	4.0675E-03	2.6	4.5231E-05	8.6	4.2066E-06	44	7.0201E-08	1657
80	8	9,345	107.3	38.0	0.423	0.039	0.00065	1.1486E-02	0.9	4.0670E-03	1.8	4.5225E-05	8.0	4.2057E-06	38	7.0050E-08	1751
90	9	9,267	106.4	37.7	0.419	0.039	0.00065	1.1487E-02	1.7	4.0676E-03	3.2	4.5233E-05	8.4	4.2071E-06	42	7.0528E-08	1725
100	10	9,182	105.5	37.4	0.415	0.039	0.00065	1.1488E-02	2.3	4.0682E-03	4.3	4.5245E-05	10.0	4.2081E-06	46	7.0671E-08	1929
								mean	1.1487E-02		4.0675E-03		4.5231E-05		4.2066E-06		7.0179E-08
								1SD	6.2081E-07		4.3349E-07		7.5843E-09		9.52215E-10		2.52131E-10
								1RSD %	54.0		106.6		167.7		226.4		3,592.7